
Ultra-Low Temperature Freezer

User Manual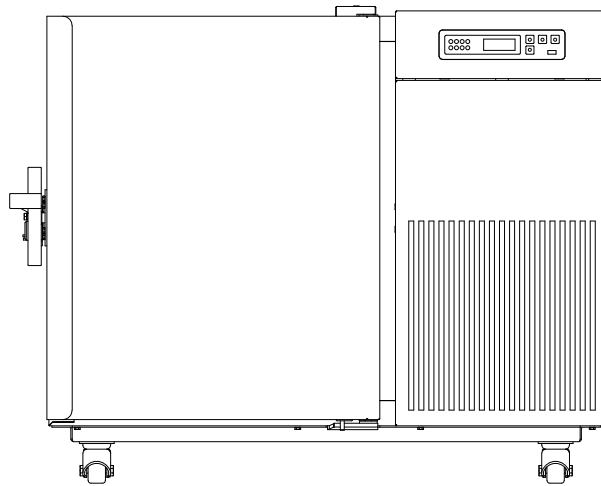

MODEL:

AULT50PS

BEFORE USE, PLEASE READ AND FOLLOW ALL SAFETY RULES AND OPERATING INSTRUCTIONS.

Models:	Serial Number:
----------------	-----------------------

Accucold Division of Felix Storch, Inc.
An ISO 9001:2015 registered company
770 Garrison Avenue
Bronx, NY 10474
www.accucold.com

Table of Contents

1. Application Notes	1
2. Safety Instructions	2
3. Precautions While in Use	4
4. Product Installation	4
4.1 Handling and Transportation	4
4.2 Installation Environment	4
4.3 Installation Site	5
4.4 Preparation Before Use	5
4.5 First Power-On	5
4.6 Operation after Power Failure	5
5. Specifications	6
6. Component Composition	7
6.1 Included with your Equipment	7
7. Operating Instructions	8
7.1 Function Description	8
7.2 Function Settings	10
7.3 Parameter Display	10
7.4 Alarm Display	11
7.5 Alarm Demonstration Function	12
7.6 Maximum and Minimum Temperature Display	12
7.7 Remote Alarm Terminal and RS485 Interface	12
8. Optional Features	13
8.1 CO ₂ Backup System	13
8.2 Freezer Rack and Freezer Box	13
9. Defrost, Discontinuation, and Maintenance	14
9.1 Equipment Maintenance	14
9.2 Equipment Discontinuation	14
9.3 Maintenance, Replacement and Recovery of Rechargeable Batteries	15
9.3.1 Battery Maintenance	15
9.3.2 Battery Replacement Procedure	15
10. Troubleshooting and Maintenance Services	16
11. Limited Warranty	21

1. Application Notes

Thank you for choosing our product. For safe, convenient operation and proper maintenance of the equipment, please read this Operation Manual carefully before use and retain it for future reference.

Operators may copy selected sections of this manual for internal use only, such as for training or emergency procedures. These sections are clearly identified in the manual's table of contents.

The manufacturer assumes no responsibility for equipment damage or malfunction resulting from failure to operate the equipment in accordance with the instructions and methods specified in this manual.

There are currently no known undesirable side effects associated with the use of the ultra-low temperature freezer when operated as instructed.

Due to continuous product improvements, the functions described in this manual may differ from those of the product you have purchased. In such cases, refer to the actual product and its installed features.

- Please carefully read the Attention and Safety Precautions in Section 2: Safety Instructions before operating the equipment.
- During transportation and use, avoid violent vibration or impact. Protect the freezer from rain. Store the unit in a clean, well-ventilated environment with humidity not exceeding 80% and free from corrosive gases.
- Maintenance and repair of the equipment must be performed only by the manufacturer or an agent authorized by the manufacturer.
- If the operator encounters a situation not addressed in this manual, contact the manufacturer or an authorized service agent for proper guidance.
- Failure to operate the equipment in accordance with the specifications and instructions in this manual may result in damage to the equipment.
- Use only accessories provided or approved by the manufacturer. The manufacturer assumes no responsibility for adverse consequences resulting from the use of unapproved accessories.
- Inspect and maintain the equipment regularly to ensure proper and reliable operation.
- The Ultra-Low Temperature Freezer is intended for cryopreservation of plasma, cells, tissues, and similar materials. It is suitable for use in blood centers, hospitals, epidemic prevention centers, research institutes, and similar facilities.
- Do not use mechanical devices or other means to accelerate the defrosting process, except those recommended by the manufacturer.
- Any serious incident involving the device must be reported to the manufacturer and to the applicable regulatory authority in the country of use.

Tips

Warning:

After each use, be sure to dry the water stains around the sealing strip to prevent freezing. Wear protective gloves when storing items in the equipment, and beware of frostbite!

- Always use protective equipment correctly (including clothes, gloves, goggles, etc.)
- Always practice good hygiene habits
- Each personnel is obliged to be responsible for his or her own safety

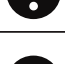
2. Safety Instructions

When using this product for the first time, please review the meanings of the following warning symbols and carefully read all safety precautions to ensure safe and proper operation of the equipment.

 Warning! Failure to observe the precautions may result in serious personal injury or death.	 Attention! Failure to observe the precautions may result in personal injury or equipment failure and related property losses
 Protective conductor terminal Must be grounded	 Risk of Explosion This sign indicates the risk of explosion when using volatile and explosive chemicals.
 Warning: Pinch point	 Warning: Low temperature
 Warning: Fire hazard	

 Warning: Failure to observe the precautions may result in serious personal injury or death.
 Do not touch the equipment with wet hands to avoid electric shock. Equipment shall not be grounded through gas pipeline, water supply pipeline, telephone line, or lightning rod, which may easily cause electric shock.
 This equipment must be installed by a qualified technician or authorized service personnel. Improper installation may result in electric shock or fire.
 Be sure to install the equipment on a solid and flat ground and take due care to prevent tipping over. If the ground is not solid enough or the installation location is not appropriate, it may cause the equipment to fall over and cause equipment damage or personal injury.
 Handle the power cord carefully to avoid short circuits or open circuits. Please turn off the power before pulling out the power plug. Hold the power plug carefully and pull it out. Do not pull the wires of the power plug. Otherwise, it may cause electric shock or fire due to the short circuit. Don't bundle the power cord, don't press it under furniture or heavy objects, and don't place it near a heat source such as the compressor.
 Insert the power plug into the outlet tightly to ensure firmness and reliability, so as to avoid electric leakage. After installation, the power plug must be within reach, so that the power plug can be unplugged in time in case of emergency.
 Separate special outlets must be used and grounded reliably. The cross-sectional area of the copper conductor in the wall connected with the outlet must be more than 4mm ² . Do not lengthen the power cord without authorization to avoid heating or fire.
 Do not use the power supply that is not specified in the equipment design, so as to avoid overheating, short circuit and other faults. For example, connecting 110V rated voltage products to 220V power supply may cause faults such as overheating and equipment burning. For detailed input voltage AC please refer to Specification (Rated voltage ±10%). If the operating voltage is too low or too high, a suitable automatic voltage stabilizer must be installed for cooperative
 Place the equipment on stable surface to avoid shaking.
 Do not place the equipment in a dangerous area, and do not operate the equipment near flammable items to prevent explosion or fire accidents.
 Vents and drains are located at the bottom of the equipment. It is forbidden to place items at the bottom of the cabinet.

	Do not place the equipment in areas exposed to the sun or rain, so as to prevent danger such as short circuit or overheating.
	Do not tilt or lay the equipment sideways, and do not impact the equipment body. Refrigeration systems are installed in the equipment, which is easy to be damaged by tilt or impact.
	Please place the equipment in a dry and dust-free environment to avoid risks such as overheating, and short circuit.
	In case of unexpected sound, smell, smoke, etc. when the power is turned on, please unplug the power in time and contact the manufacturer or supplier.
	Please place the equipment in a dry and ventilated environment, and ensure that the equipment vents and instrument surfaces are not blocked or shielded by walls or other objects. Do not use it in a poorly ventilated environment to prevent damage caused by heat released by equipment.
	It is forbidden to disassemble and modify this equipment without authorization, so as to avoid potential safety hazards. In this case, we will not bear any responsibility for quality accidents.
	It is forbidden to put inflammable and explosive dangerous goods, strong corrosive acids, alkalis and other items unsuitable for the equipment in the equipment.
	When storing toxic, hazardous, or radioactive materials, operate the equipment only in designated safe areas. Improper use may pose risks to human health and the environment.
	Metal objects such as nails or iron wires shall not be inserted into any aperture and gap or any outlet of the equipment, otherwise electric shock or injury may be caused due to accidental contact between the above objects and moving parts.
	In order to ensure the normal operation and ventilation and heat dissipation of the equipment, the back, left and right sides of the freezer body should be at least 30cm away from the wall, and the air inlet and air outlet must not.
	This equipment must be connected to a ground wire, and the equipment conductor terminal must be protected.

	Note: Failure to observe the precautions may result in personal injury or equipment failure and related property losses.
	It is forbidden to store living animals, flowers or other items with strict temperature requirements in the equipment.
	When the equipment is running, do not touch the inner surface of the cabinet without wearing protective gear.
	Hold the handle and close the door to avoid pinching your fingers. When the equipment is not used for a long time, unplug it and pack it for storage.
	When restarting the equipment after power failure or power off, check the equipment settings first, otherwise the stored items may be damaged due to the change of settings.
	The equipment can be used for item preservation, not as production equipment!
	Keep the keys in a secure location to prevent children from accidentally opening the door and causing injury.
	When handling the equipment, please be careful not to tip over the equipment, so as to prevent equipment damage or personal injury.
	When handling, it shall be lifted from the bottom, with the inclined plane not be greater than 45°, and it shall be handled with care. Use the equipment in safe areas. Improper use may cause harm to human health or environment.

3. Precautions While in Use

- When the equipment is running, the contact part between the front part of the equipment and the outer door may generate heat. This is not a fault: to prevent condensation around the casing, a heating anti-condensation pipe is installed in the equipment.
- Before placing items into the equipment, ensure that the temperature in the freezer has reached the set value first, and then put the articles in batches. Do not load more than one-third (1/3) of the freezer's internal volume at a time, as excessive loading may cause a temporary temperature rise.
- The freezer cabinet is equipped with an access port for routing a test line during testing. After the test line has been routed, the access port must be resealed with thermal insulation material. Failure to reseat the access port may prevent the freezer from reaching the set temperature and may cause condensation to form around the exterior of the access port.
- The temperature displayed on the control panel represents the temperature measured by the sensor inside the freezer. When the unit is first started, there may be a difference between the displayed temperature and the actual temperature at the center of the chamber. As the freezer reaches stable operation, the displayed temperature will gradually align with the actual internal temperature.
- Clean the equipment using a diluted neutral cleaner only. Do not use brushes, acidic cleaners, gasoline, soap powder, polishing agents, or hot water, as these may damage the painted surfaces and plastic or rubber components. Do not wipe plastic or rubber parts with volatile solvents such as gasoline.
- After a period of operation, frost may accumulate on the inner walls and inner door of the freezer. Excessive frost buildup can reduce insulation performance and increase power consumption. When the frost layer reaches approximately 5 mm, use the supplied defrosting tool to remove the frost.
- Before defrosting, remove all stored items and place them in a suitable storage environment to prevent damage caused by temperature rise during the defrosting process.
- Cooling coils are located behind and along the sides of the inner walls. Do not use sharp objects such as knives, ice picks, or screwdrivers to remove frost. Avoid scratching the inner walls during defrosting, as this may result in equipment damage or malfunction.
- If the equipment will not be used for an extended period, disconnect the power supply.

4. Product Installation

4.1 Handling and Transportation

- The product is heavy, the handling process requires the use of forklifts or hydraulic trucks inserted underneath the bottom bracket, so as not to cause damage to personnel and products.
- Avoid touching and scratching the preservation box during handling. When handling, the maximum tilt angle cannot exceed 45° (the angle between the horizontal direction), so as not to cause refrigeration system failure.
- Packaging removal: the product is packaged in a wooden bottom bracket, column packaging, and cardboard box.

4.2 Installation Environment

- Ambient temperature: 16°C ~ 32°C, the most ideal temperature is 18°C ~ 25°C, and the air conditioning system shall be used when necessary.
- Relative humidity: $\leq 80\%RH$.
- Ensure there is no strong vibration and corrosive gas around.
- Avoid areas with large amounts of dust.
- Avoid rocking or shaking the equipment.
- Elevation of the working position of the equipment: less than 2000m.
- Indoor use, pollution degree 2, and over voltage category II.
- For detailed input voltage AC refer to Specification (Rated voltage $\pm 10\%$).
- Avoid areas with direct sunlight, other cold and heat sources and strong electromagnetic interference, which will affect the normal operation of the control system and directly damage the system in severe cases.

Attention: The ambient temperature has a significant impact on equipment performance. If the specified environmental requirements are not met, the equipment may not operate properly. Ensure the environment meets the required conditions before using the equipment. The equipment is intended for intermittent operation.

4.3 Installation Site

To ensure proper operation and optimal performance, the installation site must meet the following requirements:

- Do not install the equipment in a narrow or enclosed space. The room doorway must be wide and tall enough to allow normal movement of the equipment. Adequate access must be maintained to prevent maintenance difficulties in the event of failure, which could otherwise lead to damage to stored items due to delayed repairs.
- The installation surface must be solid, level, non-combustible, and capable of supporting the full operating weight of the equipment.
- Install the equipment in a well-ventilated area and protect it from direct sunlight.
- Each unit must be connected to a dedicated power outlet. Ensure the power plug is securely connected to the outlet.
- Verify the rated operating voltage before use. In locations with unstable power supply, use an appropriate voltage stabilizer rated for the equipment load to ensure the input voltage meets installation requirements.
- The equipment must be properly grounded. If the power outlet includes a grounding connection, verify that grounding is effective before use. If the outlet is not grounded, grounding must be installed by qualified engineers or technicians before operating the equipment.

4.4 Preparation Before Use

1. Remove the outer packaging of all products (including the protective foam in the packaging box)

Warning: Don't put the plastic bags within the reach of children, so as to prevent suffocation accidents.

2. Inventory of accessories: Check the accessories and materials according to the packing list.
3. Cleaning: Clean the product once before use.
4. Adjust anti-skid leveling feet: Use a wrench to rotate the fixing foot clockwise so that the fixing foot extends outward and is supported on the ground. Ensure that the equipment will not move when in use.

4.5 First Power-on

When using the equipment for the first time, follow these steps:

1. After the equipment is placed, leveled and cleaned, it shall stand for more than 24 hours, and then power it on to ensure the normal operation of the equipment.
2. Under no-load condition, connect the power cord to a special outlet with appropriate specifications.
3. After switching on, turn on the power switch of the equipment (if the auxiliary cooling system is installed, turn off the switch first).
4. Set the required temperature in the freezer (in order to give full play to the best performance of the freezer, if the user needs to use the freezer at a storage temperature below -80°C , please set the temperature at -80°C when using it for the first time. After the system runs stably at -80°C for 3~4 hours, the temperature can be further lowered to below -80°C !).
5. In case of an alarm after power-on for the first time, click the mute button to eliminate the alarm sound.
6. Check whether the operating temperature of the equipment reaches the required value, observe the normal start and stop of the equipment for more than 24 hours, and put a small amount of items in the freezer after confirming the normal performance.
7. Please store items in batches, with the items not exceeding 1/3 of the freezer volume each time. Ensure that the equipment is running properly after shutdown for more than 12 hours before putting in the next batch of items).
8. Try not to open the door during cooling, otherwise the temperature will rise.
9. Turn on the auxiliary cooling system if it is installed.

4.6 Operation after Power Failure

1. The equipment has a memory function for the set value. When the power is restored after power failure, the equipment will continue to operate before power failure.
2. Once the equipment is powered off, it requires 5 minutes before it can be switched on again, so as to avoid damaging the compressor.
3. We guarantee the normal operation of this equipment under certain conditions, but we are not responsible for any loss or damage of stored items after power failure.

Notes:

- The equipment is designed for transportation and relocation using its integrated casters.
- A designated person should be responsible for daily monitoring and recording of the equipment's operating status (recommended inspection every 2–4 hours). In the event of a malfunction or shutdown, the internal temperature will rise. If the unit cannot be repaired promptly, remove the stored items and transfer them to a storage environment that meets their required temperature conditions to prevent damage.
- Before placing items into the equipment, confirm that the operating temperature range of the unit meets the temperature requirements of the stored materials. Failure to do so may result in damage to stored items. When loading the unit, ensure that the air inlet and air outlet are not blocked.
- Due to refrigeration system inertia, a difference may exist between the displayed temperature and the set temperature. This is a normal operating condition.
- This equipment is intended for storage only and must not be used for routine production operations. Do not load large quantities of warm items at one time. Excessive thermal loading may cause prolonged compressor operation and overheating, potentially resulting in compressor damage. Items must be loaded in batches to allow the freezer to cool gradually until the required storage temperature is reached.
- Do not use electrical appliances inside the equipment unless they are approved and certified for such use.
- During operation, short-term power supply fluctuations or brief alarm conditions will not affect normal freezer operation if the voltage variation remains within $\pm 10\%$ of the rated value.
- Do not change the set temperature frequently within short time intervals, as temperature inertia may prevent the system from achieving the desired setpoint. When loading items, ensure adequate air circulation space around the cabinet interior, and do not block the internal temperature sensor, as this will affect temperature stability and control accuracy.
- When loading items, ensure they are properly sealed. Excessive or insufficient moisture content in stored items may affect humidity levels inside the cabinet. Ambient humidity and environmental conditions also influence internal humidity, particularly if the door is opened frequently or not fully closed.


Warning: Children must not play with this equipment. Failure to comply may result in injury or damage.

5. Specifications

Model	Ambient temperature	Climate type	Refrigerant	(Rated) voltage (v~)	Rated frequency (Hz)	Temp Range (°C)	Volume (L)	(Rated) current (A)	Net Weight (kg)	Exterior dimensions (DxWxH) (mm)
AULT50PS	16~32°C	N	Mixture Gas	110	60	-40~-86	50	10	96	700X951X757

* The foaming material of this product is cyclopentane. If optional electromagnetic lock module is selected the height of the whole unit is increased by 45mm.

6. Component Composition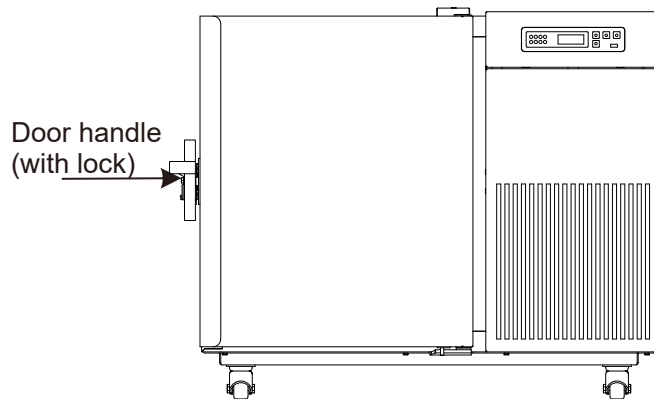


* Due to ongoing product improvements and model variations, the actual product may differ from the illustrations shown. Refer to the actual product for details. Illustrations are provided for functional reference only.

Structure and Composition:

The product consists of a freezer cabinet, door (glass door or insulated foam door), refrigeration system, and control system.

Opening the Door: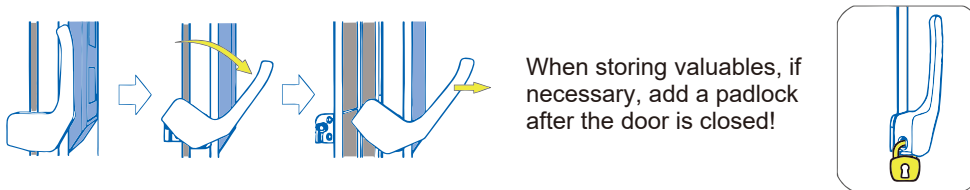

Note: For products equipped with an electric control lock safety function, log in to the system and release the electric control lock before attempting to open the door.



6.1 Included with your Equipment

Name	Operation Manual	Certificate of Conformity	USB flash drive	Key	Defrosting shovel
Number	1	1	1	2	1

* The specific accessories shall be subject to the physical objects received.

7. Operating Instructions

7.1 Function Description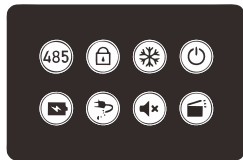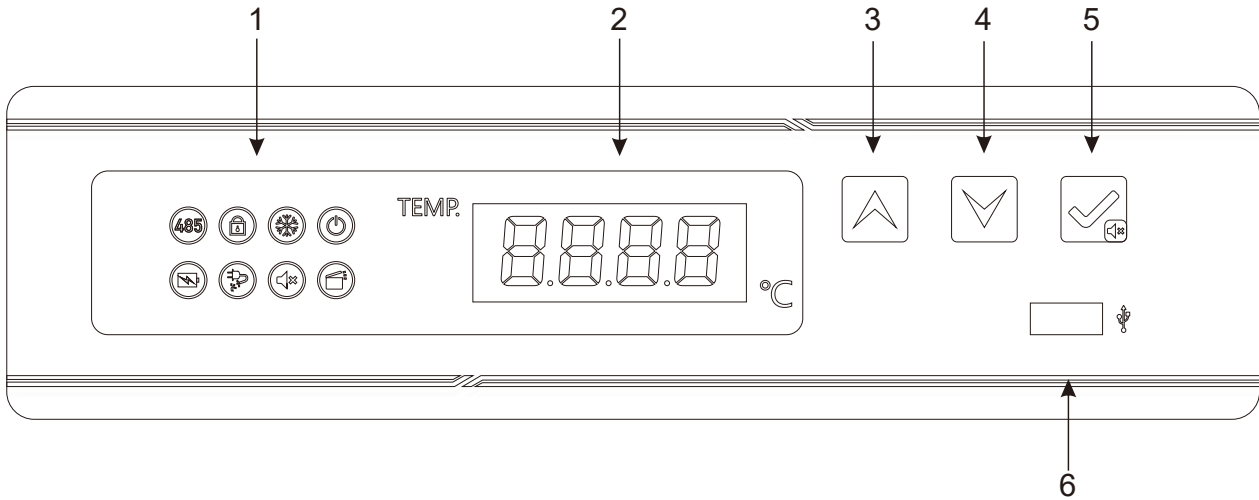

a. Remote Port	b. Key Lock	c. Refrigeration Indicator	d. Power Indicator
e. Low Battery Level Indicator	f. Power Failure	g. Mute Indicator	h. Door Switch Indicator

1. Function Indicators (see table above)

a. Remote Port Indicator

- When the reserved RS-485 remote port is not connected, the remote port indicator remains off.
- When the freezer is successfully connected to the RS-485 serial port, the indicator turns on.

b. Key Lock Indicator

When the keys are locked, none of the buttons will respond and the key lock indicator will illuminate. To unlock the keys:

1. Press and hold the **Up** and **Down** keys simultaneously for 3 seconds.
2. Enter the password (default: **0005**).
3. Press the **Set/Mute** key to confirm.
4. The key lock will be released and the indicator will turn off.

In the unlocked state, if no key is pressed for 60 seconds, the key lock will activate automatically, and the indicator will turn on.

To manually lock the keys, press and hold the **Up** and **Down** keys for 3 seconds.

c. Refrigeration Indicator

- If the compressor is in working condition, the refrigeration indicator is normally on.
- If the compressor is in a shutdown state, the refrigeration indicator is off.

d. Power Indicator

When the indicator is on, it means that the power switch of the product has been turned on.

e. Low Battery Level Indicator

- When the battery voltage is less than 10.8V, the buzzer will sound, the low battery indicator will turn on, and the low battery code **[BL]** will flash on the screen.
- When the battery voltage is greater than 12V, the buzzer will stop, the low battery indicator will turn off, and the screen will return to normal display.

f. Power Failure

The freezer is normally powered by 220V. When there is a power outage, the buzzer will sound, the screen will flash the power failure code **[PF]**, and the power-off indicator will turn on. When the input power is turned on, it will return to normal and the power-off indicator will turn off.

g. Mute Indicator

- When the mute key is activated, the indicator is on.
- When the mute function is canceled, the indicator is off.

h. Door Switch Indicator

- When the door is opened, the door switch indicator is on.
- When the door is closed, the door switch indicator is off.

2. Temperature Display Window:

The temperature display window/screen will display the average temperature inside the freezer compartment in Celsius (°C) under normal operations.

View ambient temperature:

- During normal operation, press **Set/Mute** and the screen will display the ambient temperature, and return to normal display after 5 seconds of inactivity.
- In the key locked state, press **Set/Mute** and **Down** and the screen will display the ambient temperature. The screen will return to normal after 5 seconds of inactivity or press **Up** and **Down** simultaneously.

3. Up:

In parameter setting, press **Up** to move to the next parameter or increase the parameter value.

- When adjusting the set temperature, press **Up** to increase the set temperature value.
- When adjusting the parameter value, press and hold **Up** and the parameter will increase rapidly.

During normal operation, press and hold **Up** to import 12 months of USB data.

4. Down:

In parameter setting, press **Down** to move to the previous parameter or decrease the parameter value.

- When adjusting the set temperature, press **Down** to reduce the set temperature.
- When setting the parameter value, press and hold **Down**, and the parameter will decrease rapidly.

5. Set/Mute:

During normal operation press **Set/Mute** and display the ambient temperature for 5 seconds. The display will then return to the freezer temperature and alarm status.

Press and hold **Set/Mute** to enter the user menu.

When the alarm is activated and keys are unlocked (examples: high-temperature alarm, door opening alarm, sensor failure alarm, etc.) press **Set/Mute** to silence the alarm and the ambient temperature will display for 5 seconds, then the return to the freezer compartment temperature and alarm status.

Note: Muting only silences the alarm for the current alarm condition. If another abnormal condition occurs, the alarm will sound again.

Press the **Set/Mute** again to reactivate the alarm, the ambient temperature will display for 5 seconds and then return to freezer compartment temperature and alarm status.

During normal operation (keys unlock), **Set/Mute** can be used as a setting key.

In parameter setting and keys are unlocked, press **Set/Mute** to display parameter names and values. Press and hold **Set/Mute** to save the settings and return to the normal display.

6. USB Data Export:

Automatic Export: When a USB flash drive is inserted into the USB port, the unit will beep once and the screen will display **[on]**. PDF files containing data that has not yet been exported will be automatically generated and saved to the USB drive. After data transfer is completed, the unit will beep again and the screen will display **[End]**. After 6 seconds, it will return to normal interface.

Note: If the amount of data is small, the display may not show **[on]** and **[End]** during the export process.

Manual Export: During normal operation, when the USB flash drive is connected and the file is not being generated, press and hold **Up** and the screen will display **[d01]**.

Press **Up** or **Down** to select:

[d00] – Generate the current data file.

[d01–d12] – Generate the PDF file for record data from previous months (1–12).

Press the **Set/Mute** to confirm and begin file generation.

Note: When the alarm of the screen flashes **[LoF]** the recorder is not started. To activate, press and hold **Set/Mute** and **Up** simultaneously till **[LoF]** disappears. A beep will sound to confirm that the recorder has started.

7.2 Function Settings

1. Power On

Power on the unit and the power switch, the equipment will enter the working state.

2. User Parameter Settings:

To unlock parameter settings under normal operation:

1. Press and hold **Up** and **Down** simultaneously and the screen will display the parameter code [0000].
2. Press **Up** to enter the password [0005].
3. Press and hold **Set/Mute** to confirm and unlock.
4. Press and hold **Set/Mute** again, the screen will display the parameter code [PS1] and enter the parameter settings.
5. Use the **Up** or **Down** key to scroll the parameters.

To adjust parameter settings:

1. Press **Set/Mute** to enter the parameter group, and the first parameter name in that group will be displayed.
2. Use **Up** or **Down** to scroll the parameters.
3. Press **Set/Mute** to display the current value of the selected parameter.
4. Use **Up** or **Down** to adjust the value.
5. Press **Set/Mute** to temporarily store the new value and return to the parameter list.
6. Repeat steps 1 – 5 to adjust additional parameters as needed.
7. Press and hold the **Set/Mute** key to save the modified parameters and return to the parameter group menu.
8. Press and hold **Set/Mute** again or do not press any key for 1 minute to exit the parameter setting program.

7.3 Parameter Display

The following table is a comparison table for parameter display.

No.	Menu item	Parameter Range	Suggested settings	Remarks
1	MAX	—	—	The highest temperature since last clearance
2	MIN	—	—	The lowest temperature since last clearance
3	CLR	—	—	Clearance of the Max and Min temperature records
4	CF	C/F	—	Temperature unit selection, C° Celsius, F° Fahrenheit
4	Set	-86.0~-40.0	-80.0	Temperature setting
5	H	0.0~10.0	10.0	Set value of high temperature alarm set+H; When H =0, High temp alarm is disabled; When the alarm is over high temp alarm set, H1 will be displayed on the controller
6	L	0.0~10.0	10.0	Set value of low temperature alarm set-L; When L =0, Low temp alarm is disabled; When the alarm is below low temp alarm set, L1 will be displayed on the controller
7	n	Set logger module time - year	—	—
8	y	Set logger module time - month	—	—
9	r	Set logger module time - day	—	—
10	s	Set logger module time - hour	—	—
11	F	Set logger module time - minute	—	—
12	Pt	0~240min	20	Print interval
13	tH1	20.0~50.0°C	40.0	Upper limit of ambient temperature alarm
14	Ps1	0~9999	5	User menu password settings
15	b1	—	—	Repair Information 1
16	b2	—	—	Repair Information 2

Quick setting of time after power on

After the power-on self-test on the display board is completed, the quick setting menu is displayed.

Menu item	Menu	Menu description	Set range	Default	Unit
Quick Settings menu	n	Set logger module time - year	10~50		/
	y	Set logger module time - month	1~12	--	/
	r	Set logger module time - day	01~31	--	/
	S	Set logger module time - hour	00~23	--	/
	F	Set logger module time - minute	00~59	--	/
	Pt	Print interval	0~240	20	min
	SCY	Temperature data recording period	0~240 0: shielded recorder	10	min

If there is no operation for 60 seconds under the quick setting menu, it will automatically exit the quick setting menu and return to normal display.

7.4 Alarm Display

The following table is a comparison table of alarm codes and fault descriptions.

Alarm Code	Default Description
H1	High temperature alarm
L1	Low temperature alarm
H2	High ambient temperature alarm
H3	High condenser temperature alarm
do	Door opening alarm
PF	Power failure alarm
bL	Low battery alarm
ER	The recorder is not connected
LoF	The recorder is not started
EE	Communication failure

7.5 Alarm Demonstration Function

1. In the key lock state, press and hold **Set/Mute**.
2. The display will show the temperature value and alarm code for the over-temperature alarm.
3. The alarm will sound.

7.6 Maximum and Minimum Temperature Display

When the keys are locked:

Press Up to view the maximum recorded temperature.
After blinking three times, the maximum value will be displayed.

Press Down to view the minimum recorded temperature. After blinking three times, the minimum value will be displayed.

To clear the recorded maximum and minimum values:


Press and hold Up + Down + Set/Mute at the same time.
The buzzer will beep to confirm the data has been cleared.

! Notes:

- It is normal for a High Temperature Alarm to activate when the unit is first powered on. Load items into the freezer only after the temperature has stabilized and the alarm clears automatically.
- Do not load more than 1/3 of the freezer capacity at one time. After the unit reaches the set temperature and operates normally for 24 hours, you may load an additional 1/3 of the capacity.
- Avoid opening the door during the cooling process, as this will cause the internal temperature to rise.
- Do not change the set temperature frequently within a short period. Due to temperature inertia, frequent adjustments may prevent the unit from reaching the desired temperature.
- Ensure there is adequate air circulation inside the cabinet when loading items. Do not block the temperature sensor opening, as this may affect temperature stability and accurate display.

7.7 Remote Alarm Terminal and RS485 Interface

The remote alarm terminal is installed on the back or the lower part of the right side of the freezer body, and the alarm signal is output by the terminal. The load-bearing capacity of the terminal is 30V, 2A DC. Users can choose normally open (N.O.), normally closed (N.C.) and common terminals (COM) according to their own needs.

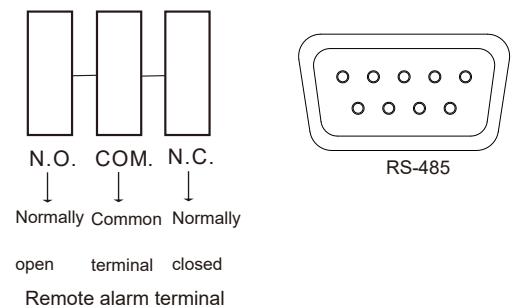8. Optional Features

8.1 CO₂ Backup System

The CO₂ backup system is an auxiliary refrigeration system that uses liquid CO₂ for refrigeration. It provides refrigeration assistance for the equipment in case of power failure or refrigeration failure, to ensure that the temperature in the equipment is maintained within the required temperature range for a certain period of time.

When the equipment works normally, the CO₂ backup system is in standby state. When the equipment is powered off, the system uses its own rechargeable battery to work and spray liquid CO₂ into the refrigeration equipment at intervals to achieve the refrigeration effect. In case of refrigeration failure of the equipment, the forced operation switch can control the backup system to work forcibly to provide refrigeration.

Notes:

- Ensure the cylinder is filled with liquid carbon dioxide (CO₂) .
- Liquid CO₂ is extremely cold. Use caution during installation and operation to prevent frostbite.
- The CO₂ backup system is intended for emergency use only. Test the system once per month to ensure proper operation.
- Install and operate the unit in a well-ventilated area to prevent discomfort caused by excessive carbon dioxide or reduced oxygen levels.

8.2 Freezer Rack and Freezer Box

Model	Internal dimensions of the freezer body W x D x H	Shelf/inner door	Number of freeze racks	Dimensions of the freeze rack W x D x H	Number of freeze boxes that can be placed (2 in.)
AULT50PS	12" x 16 ¾" x 16 ⅞" (305 x 425 x 430)	1/1	2 pcs	5 ½" x 16 ¼" x 13 ⅞" 140 x 413 x 351	36 pcs

NOTE: For assistance with optional features, contact us at **1-888-4-MEDLAB**, Monday through Friday, 9:00 AM–5:00 PM ET, or visit our website at **www.accucold.com/support** anytime for ordering and installation information.

To order replacement parts, visit our website: **summitapplianceparts.com**

9. Defrost, Discontinuation, and Maintenance

9.1 Equipment Maintenance

Defrosting

When frost buildup on the inner walls becomes excessive, cooling performance may decrease and power consumption may increase.

- Use the supplied defrosting shovel to remove frost.
- Do not use sharp or metal tools, as they may damage the unit.
- Before defrosting, disconnect the power supply and remove all stored items.
- After defrosting, restore power and allow the unit to reach the set temperature before placing items back inside.
- Always wear protective gloves when removing or loading items to prevent frostbite.

Cleaning and General Maintenance

- Clean the unit regularly.
- Before cleaning, unplug the power cord for safety.
- Wipe the interior and exterior surfaces with a soft cloth dampened with warm water.

Cleaning the Filter Screen

Regularly remove the dust on the filter screen to ensure the desirable heat dissipation effect.

Warning:

- Disconnect the power supply before performing any repair or maintenance to prevent electric shock or injury.
- Avoid inhaling any harmful substances or airborne particles during maintenance, as this may pose health risks.

Notes:

- Do not spray water directly onto the cabinet, as this may reduce electrical insulation and cause metal parts to rust.
- Do not use hot water, corrosive detergents, or organic solvents to clean the unit.
- Do not place heavy objects on top of the equipment, as this may cause deformation.
- The built-in air switch supports both manual operation and automatic protection. In the event of a short circuit or overload, it will automatically disconnect the power supply to protect the unit and electrical system.

9.2 Equipment Discontinuation

Vacation

If the equipment will not be used for an extended period or is stored in an unsupervised area:

- Disconnect the power supply.
- Clean the interior and exterior surfaces with a soft cloth dampened with warm water.
- Allow the unit to dry completely.
- Close and seal the unit properly.
- Lock the equipment to prevent children from opening the door.


Disposal

When the equipment reaches the end of its service life, it must be disposed of in accordance with local regulations. The unit should be delivered to a qualified recycling or disposal agency. Do not allow unqualified personnel to dismantle or disassemble the equipment. Discarded equipment must be placed in a designated area that is inaccessible to children to prevent injury.

9.3 Maintenance, Replacement and Recovery of Rechargeable Batteries

9.3.1 Battery Maintenance

1. To extend battery life, operate the unit for at least 24 hours each month to ensure proper charging.
2. If the equipment is not used for an extended period, connect it to the main power supply at least once a month. Turn on the power switch and allow the unit to run for no less than 24 hours to fully charge the battery.
3. If the main power supply is interrupted, turn off the power switch immediately to prevent battery discharge, which may cause permanent damage.
4. During long-term power outages or transportation, ensure the main power switch is turned off. Prolonged discharge may result in battery damage and abnormal display when power is restored.
5. The battery is a consumable component with an expected service life of approximately 2 – 3 years. Improper use or end-of-life conditions may trigger a low battery alarm. Refrigeration performance will not be affected; however, alarm and power-failure functions may not operate properly. Contact authorized service personnel for battery replacement if needed. For assistance call our Customer Service facility at **800-932-4267** (Ext. **513**) between 9:00 AM and 5:00 PM ET or visit our website **www.accucold.com/support** at any time.

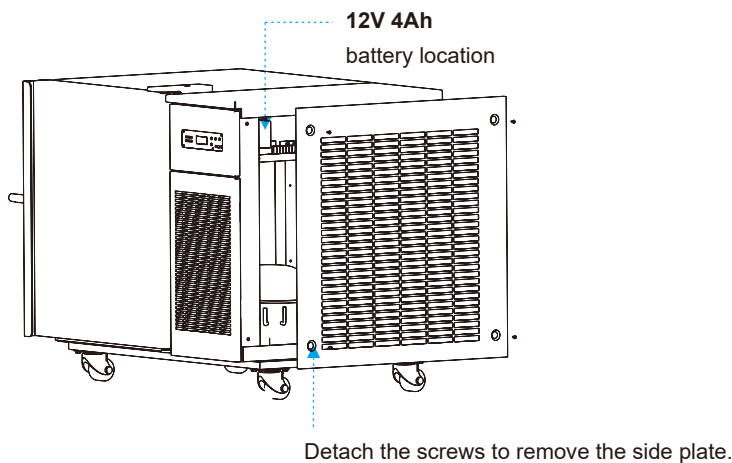9.3.2 Battery Replacement Procedure

1. Turn off the power switch and unplug the unit from the electrical outlet.
2. Using a screwdriver, remove the screws on the right-side panel of the freezer frame.

Note:

The battery can be replaced without opening the electrical box. The electrical box contains high-voltage components. If it must be opened, disconnect the power supply, unplug the unit, and turn off the power switch. The electrical box may only be opened by qualified technicians or authorized service personnel.

3. Locate the battery on the left side outside the electrical box and disconnect the battery cables.

- Before disconnecting, note the polarity and cable connections.
- Incorrect connection of the positive and negative terminals may damage the control system.
- Typically, the red wire connects to the positive terminal and the black wire connects to the negative terminal.

4. Remove the screws securing the battery fixing plate and carefully remove the battery.
5. Replace with a battery of the same specification: **BT-12M4.0 AC (12V, 4.0Ah)**.
6. Dispose of the used battery in accordance with local regulations. Contact a local battery recycling agency for proper disposal.

To ensure correct installation and prevent damage to the control system, contact authorized after-sales service personnel for battery replacement or guidance.

10. Troubleshooting and Maintenance Services

You can solve many common problems easily, saving you the cost of a possible service call. Try the suggestions below to see if you can solve the problem before calling the servicer.

Problems	Causes and Solutions
Ultra-low temperature freezer does not operate	<ul style="list-style-type: none"> • Ensure the outlet is powered. • Make sure the power plug is securely connected. • Check that the fuse is not blown or disconnected. • Verify that the supply voltage is within the specified range (not too high or too low).
Compressor is not running	<ul style="list-style-type: none"> • Ensure the power switch (located on the lower right side of the unit) is turned on. • Verify that the temperature setting is correct.
The temperature does not reach the set value	<ul style="list-style-type: none"> • Check that frost buildup on the inner walls is not excessive. • Ensure the door is fully closed and avoid frequent door openings. • Do not overload the freezer at one time. • Make sure the ambient temperature is not too high. • Check that the filter screen is clean and not blocked.
High noise	<ul style="list-style-type: none"> • Ensure the unit is placed on a flat and stable surface. • Make sure the unit is not touching the wall or other objects.
High ambient temperature alarm	<p>Cause: Ambient temperature exceeds 89.6°F (32°C) or the unit is placed too close to the wall, restricting ventilation.</p> <p>Solution: Maintain at least 12 inch (30 cm) clearance between the unit and the wall. Keep ventilation openings clear. Clean the filter screen regularly. Improve room ventilation or use air conditioning during hot weather.</p>

The Following Conditions Are Normal

1. A slight clicking sound when the compressor starts or stops.
2. The sound of flowing refrigerant, similar to running water.
3. Light frost on the cabinet surface during humid or rainy weather. This can be wiped off with a dry cloth.
4. If the door is opened again within 5 minutes, it may be difficult to reopen due to internal pressure differences.
5. Minor cracks around the inner liner opening caused by temperature variations. This is normal and does not affect operation.
6. The LCD display may flash during printing or briefly during startup while data is being processed. This is normal. If the display does not return to normal, turn off the power and restart the unit.

If you have checked the information above and you still need help with your appliance, call our Customer Service facility at **1-888-4-MEDLAB** between 9:00 AM and 5:00 PM ET or visit our website **www.accucold.com/support** at any time. We will do our best to answer your questions.

To order replacement parts, visit our website: **summitapplianceparts.com**

Warning:

This product uses a flammable refrigerant. Do not disassemble or repair the unit without authorization, especially the compressor, refrigeration system, or related components. In the event of a malfunction, contact our authorized Service Center for repair.

Notes:

- Repairs, maintenance, or modifications to the equipment must be performed only by engineers certified by our company to ensure proper operation and compliance with applicable safety standards.
- Before contacting a maintenance engineer, clean and disinfect the equipment. During the warranty period, the Company will not be responsible for faults or damage caused by improper use.
- **Storage Conditions:**
Ambient temperature: -40°C to +55°C
Relative humidity: 10% to 90%

WARNING: This product can expose you to chemicals including Nickel (Metallic) which is known to the State of California to cause cancer.

For more information go to www.P65Warnings.ca.gov

Note: Nickel is a component in all stainless steel and some other metallic compositions.

CALIFORNIA CARB/SNAP DISCLOSURE

This product uses eco-friendly hydrocarbon refrigerant and fully complies with California CARB regulations.

However, we are required by California Law to provide the following disclosure statement in every product sold in California.

"This equipment is prohibited from use in California with any refrigerants on the 'List of Prohibited Substances' for that specific end-use, in accordance with California Code of Regulations, title 17, section 95374. This disclosure statement has been reviewed and approved by Felix Storch, Inc. and Felix Storch, Inc. attests, under penalty of perjury, that these statements are true and accurate."

This product does **not** use any refrigerants on the 'List of Prohibited Substances'"

NOTES

NOTES

NOTES

LIMITED WARRANTY

FIVE-YEAR LIMITED WARRANTY

Warranty coverage is available within the 48 contiguous United States, along with select areas of Alaska, Hawaii, and Canada. For locations not covered, consumers can either contact a local provider or coordinate with Felix Storch, Inc. to find nearby service options, which may qualify for partial reimbursement.

The warranty for this appliance is valid for five years from the date of purchase, provided the appliance is operated and maintained according to the instructions included with the product. The warrantor will cover the cost of factory-specified parts and repair labor needed to correct defects in materials or workmanship. Service must be performed by a designated service company.

All parts, except the compressor, are warranted for five years against manufacturing defects. Plastic parts, shelves, and cabinets are produced to commercially acceptable standards but are not covered for damages incurred during handling or for breakage.

12-YEAR COMPRESSOR WARRANTY

1. The compressor is covered for 12 years.
2. Replacement does not include labor.

ITEMS WARRANTOR WILL NOT PAY FOR:

1. Service calls to correct the installation of your appliance, to instruct you how to use your appliance, to replace or repair fuses or to correct wiring or plumbing.
2. Service calls to repair or replace appliance light bulbs or broken shelves. Consumable parts (such as filters) are excluded from warranty coverage.
3. Damage resulting from accident, alteration, misuse, abuse, fire, flood, acts of God, improper installation, installation not in accordance with electrical or plumbing codes, or use of products not approved by warrantor.
4. Replacement parts or repair labor costs for units operated outside the 48 contiguous United States, along with select areas of Alaska, Hawaii, and Canada.
5. Repairs to parts or systems resulting from unauthorized modifications made to the appliance.
6. Expenses for travel and transportation for product service in remote locations.
7. The removal and reinstallation of your appliance if it is installed in an inaccessible location or is not installed in accordance with published installation instructions.

DISCLAIMER OF IMPLIED WARRANTIES; LIMITATION OF REMEDIES

CUSTOMER'S SOLE AND EXCLUSIVE REMEDY UNDER THIS LIMITED WARRANTY SHALL BE PRODUCT REPAIR AS PROVIDED HEREIN. IMPLIED WARRANTIES, INCLUDING WARRANTIES OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE, ARE LIMITED TO FIVE YEARS. WARRANTOR SHALL NOT BE LIABLE FOR INCIDENTAL OR CONSEQUENTIAL DAMAGES. SOME STATES DO NOT ALLOW THE EXCLUSION OR LIMITATION OF INCIDENTAL OR CONSEQUENTIAL DAMAGES, OR LIMITATIONS ON THE DURATION OF IMPLIED WARRANTIES OF MERCHANTABILITY OR FITNESS, SO THESE EXCLUSIONS OR LIMITATIONS MAY NOT APPLY TO YOU. THIS WARRANTY GIVES YOU SPECIFIC LEGAL RIGHTS AND YOU MAY ALSO HAVE OTHER RIGHTS, WHICH VARY FROM STATE TO STATE.

accucold®

Accucold Division of Felix Storch, Inc.

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