accucold®

 **BOD** REFRIGERATED INCUBATOR
PERFORMANCE CHAMBER

BOD Refrigerated Incubator

ABOD6

ABOD15

ABOD18

User Manual

BEFORE USE, PLEASE READ AND FOLLOW ALL SAFETY RULES AND OPERATING INSTRUCTIONS

Write Model and Serial Numbers here:

Model _____

Serial No. _____

Accucold Division of Felix Storch, Inc.
An ISO 9001:2015 registered company
770 Garrison Avenue
Bronx, NY 10474
www.accucold.com

EQUIPMENT SAFETY

Your safety and the safety of others are very important.

We have provided many important safety messages in this manual and on your unit. Always read and obey all safety messages.

This is the Safety Alert Symbol. The symbol alerts you to potential hazards that can kill or injure you and others. All safety messages will follow the Safety Alert Symbol and either the words "DANGER" or "WARNING".

DANGER means that failure to heed this safety statement may result in severe personal injury or death.

WARNING means that failure to heed this safety statement may result in extensive product damage, serious personal injury, or death.

All safety messages will alert you about the potential hazard, tell you how to reduce the chance of injury, and let you know what can happen if the instructions are not followed.

EQUIPMENT SAFETY

Before the equipment is used, it must be properly positioned and installed as described in this manual, so read the manual carefully. To reduce the risk of fire, electrical shock or injury when using this equipment, follow basic precautions, including the following:

- Plug into a grounded 3-prong outlet, do not remove grounding prong, do not use an adapter, and do not use an extension cord.
- Replace all panels before operating.
- It is recommended that a separate circuit serving only your unit be provided. Use receptacles that cannot be turned off by a switch or pull chain.
- Never clean the equipment parts with flammable fluids. These fumes can create a fire hazard or explosion. And do not store or use gasoline or other flammable vapors and liquids in the vicinity of this or any other pieces of equipment. The fumes can create a fire hazard or explosion.
- Before proceeding with cleaning and maintenance operations, make sure the power line of the unit is disconnected.
- Do not connect or disconnect the electric plug when your hands are wet.
- Unplug the unit or disconnect power before cleaning or servicing. Failure to do so can result in electrical shock or death.
- Do not attempt to repair or replace any part of your unit unless it is specifically recommended in this manual. All other servicing should be referred to a qualified technician.
- These units are CFC- and HFC-free and contains a small quantity of Isobutane (R600a), and the ABOD18 are CFC- and HFC-free and contain a small quantity of propane (R290). Both are environmentally friendly, but flammable. They do not damage the ozone layer, nor do they increase the greenhouse effect. Care must be taken during transportation and setting up of a unit that no parts of the cooling system are damaged. Leaking coolant can ignite and may damage the eyes.
- In the event of any damage:
 - Avoid open flames and anything that creates a spark,
 - Disconnect from the electrical power line,
 - Air the room in which the unit isolated for several minutes, and
 - Contact the Service Department for advice.

- The more coolant there is in a unit, the larger the room it should be installed in. In the event of a leakage, if the unit is in a small room, there is the danger of combustible gases building up. For every ounce of coolant, at least 325 cubic feet of room space is required. The amount of coolant in the unit is stated on the data plate inside the unit. It is hazardous for anyone other than an Authorized Service Person to carry out servicing or repairs to this piece of equipment.
- Take serious care when handling, moving, and using the unit to avoid either damaging the refrigerant tubing or increasing the risk of a leak.
- Replacing component parts and servicing shall be done by factory authorized service personnel so as to minimize the risk of possible ignition due to incorrect parts or improper service.

⚠ WARNING ⚠

FOLLOW WARNING CALLOUTS BELOW ONLY WHEN APPLICABLE TO YOUR MODEL

- Use two or more people to move and install **unit**. Failure to do so can result in back or other injury.
- To ensure proper ventilation for your **unit**, the front of the unit must be completely unobstructed. Choose a well-ventilated area with temperatures above 60°F (16°C) and below 90°F (32°C). [For optimal performance, install the unit where the ambient temperature is between 72° and 78°F (23°- 26°C).] This unit must be installed in an area protected from the elements, such as wind, rain, waterspray or drips.
- The unit should not be located next to ovens, grills or other sources of high heat.
- The unit must be installed with all electrical, water and drain connections in accordance with state and local codes. A standard electrical supply (115 V AC only, 60 Hz), properly grounded in accordance with the National Electrical Code and local codes and ordinances, is required.
- Do not kink or pinch the power supply cord of the unit.
- The size of the fuse (or circuit breaker) should be 15 amperes.
- It is important that the equipment be leveled in order to work properly. You may need to make several adjustments to level it.
- All installations must be in accordance with local plumbing code requirements.
- Make certain that the pipes are not pinched, kinked or damaged during installation.
- Check for leaks after connection.
- Never allow children to operate, play with or crawl inside the unit.
- Do not use solvent-based cleaning agents or abrasives on the interior. These cleaners may damage or discolor the interior.
- Use this equipment only for its intended purpose as described in this *Instruction Manual*.
- Keep fingers out of the “pinch point” areas. Clearances between the door and cabinet are necessarily small. Be careful closing the door when children are in the area.

⚠ DANGER ⚠ Risk of child entrapment!

Child entrapment and suffocation are not problems of the past. Junked or abandoned appliances are still dangerous, even if they will “just sit in the garage a few days”.

Before discarding your old refrigerated incubator:

- *Take off the doors*
- *Leave the shelves in place so that children may not easily climb inside.*

-SAVE THESE INSTRUCTIONS-

INSTALLATION INSTRUCTIONS

Before Using your Equipment

- Remove the exterior and interior packaging.

CAUTION: After unpacking you **MUST** allow this unit to stand upright for at least 2 hours to allow the lubricant and refrigerant to drain back into the compressor and stabilize. Failure to do so may adversely affect performance and the lifetime of this unit.

- Clean the interior surface with lukewarm water using a soft cloth.

Installation of your Equipment

NOTE: Do not store or install the unit outdoors.

- Units with **-BI** in their suffix are designed to be built-in or recessed installation. The rest of the units listed in this manual are designed for freestanding use. **All models are for indoor use only.**

CAUTION: This equipment is designed for the incubation and storage of biological and laboratory samples only. Do not store beverages, perishable food items, or non-laboratory materials in this unit.

- Place the unit on a floor that is strong enough to support it when it is fully loaded. To level the unit, adjust the front leveling legs.
- For free-standing installation, allow at least **5 inches** (127mm) of space between the back, top, and sides of the unit. This allows the proper air circulation to cool the compressor and condenser for energy saving. Even for recessed or built-in installation (units with **-BI** in the model number), it is a must to keep $\frac{1}{4}$ " (6.35mm) of space on each side, and **2 inches** (51mm) on the top and at the rear. The air vent at the front of the unit must never be covered or blocked in any way.

NOTE: It is recommended that you do not install the unit near an oven, radiator or other heating source. Direct sunlight may affect the acrylic coating and heat sources may increase electrical consumption. Don't install in a location where the temperature will fall below 60°F (16°C). For best performance, do not install the unit behind a cabinet door or block the basegrille.

- Avoid locating the unit in moist areas.
- Plug the unit into an exclusive, properly grounded wall outlet. Do not under any circumstances cut or remove the third (ground) prong from the power cord. Any questions concerning power and/or grounding should be directed toward a certified electrician or an authorized service center.

Electrical Connection

Improper use of the grounded plug can result in the risk of electrical shock. If the power cord is damaged, have it replaced by a qualified electrician or an authorized service center.

This unit should be properly grounded for your safety. The power cord of this unit is equipped with a three-prong plug which mates with standard three-prong wall outlets to minimize the possibility of electrical shock.

Do not under any circumstances cut or remove the third (ground) prong from the power cord supplied. For personal safety, this equipment must be properly grounded.

This unit requires a standard 115/120 Volt AC ~ 60Hz three-prong grounded electrical outlet. Have the wall outlet and circuit checked by a qualified electrician to make sure the outlet is properly grounded. When a standard 2-prong wall outlet is encountered, it is your responsibility and obligation to have it replaced with a properly grounded 3-prong wall outlet. To prevent accidental injury, the cord should be secured behind the unit and not left exposed or dangling.

The unit should always be plugged into its own individual electrical outlet which has a voltage rating that matches the rating label on the unit. This provides the best performance and also prevents overloading house wiring circuits that could cause a fire hazard from overheating. Never unplug the unit by pulling on the power cord. Always grip the plug firmly and pull straight out from the receptacle. Repair or replace immediately all power cords that have become frayed or otherwise damaged. Do not use a cord that shows cracks or abrasion damage along its length or at either end. When moving the equipment, be careful not to damage the power cord.

Electrical Connection

Because of potential safety hazards under certain conditions, it is strongly recommended that you do not use an extension cord with this equipment. However, if you must use an extension cord it is absolutely necessary that it be a UL/CUL-Listed, 3-wire grounding type equipment extension cord having a grounding type plug and outlet and that the electrical rating of the cord be 115 volts and at least 10 amperes.

Reversing the Door

Unless ordered as the **LHD** option, your unit comes with a right-hand self-closing door swing by default. If you would like to reverse the door to open in the opposite direction, please follow the steps below:

NOTE: Reversing the door will undo the self-closing function. Doors on units with -DL2B data logger suffix are not reversible.


1. Carefully lay the unit on its back and remove the two (2) front leveling legs
2. Remove the two (2) hex head screws that connect the door hinges to the body on the bottom of the unit
3. Remove the two (2) flathead screws that connect the hinge assembly to the bottom of the door
4. Remove the white hinge from the square rod
5. Slide the door towards the bottom until it is free of the top hinge
6. Unscrew the top hinge pin from the unit and reinstall on the left side
7. Flip the door so the handle is now on the right side
8. Push the door onto the top hinge piece
9. Reinstall the hinge assembly on the left side bottom of the door using the flathead screws
10. Push the white hinge onto the square rod at a position of 9 o'clock
11. Turn the white hinge clockwise until the holes line up with screws located on the bottom of the unit
12. Replace the hex head screws and front leveling legs

NOTE: Once the door has been reversed, the BOD Refrigerated Incubator sticker will be upside down. You can remove and reapply to the preferred position. If you would like a replacement sticker, contact our parts department at parts@summitappliance.com

CAUTION: After positioning the unit, you **MUST** allow it to stand upright for at least 2 hours to allow the lubricant and refrigerant to drain back into the compressor and stabilize. Failure to do so may adversely affect performance and the lifetime of this unit.

Stabilization Period - Prior to use, the unit shall be allowed to run and stabilize for at least twenty-four hours.

LOCATION OF PARTS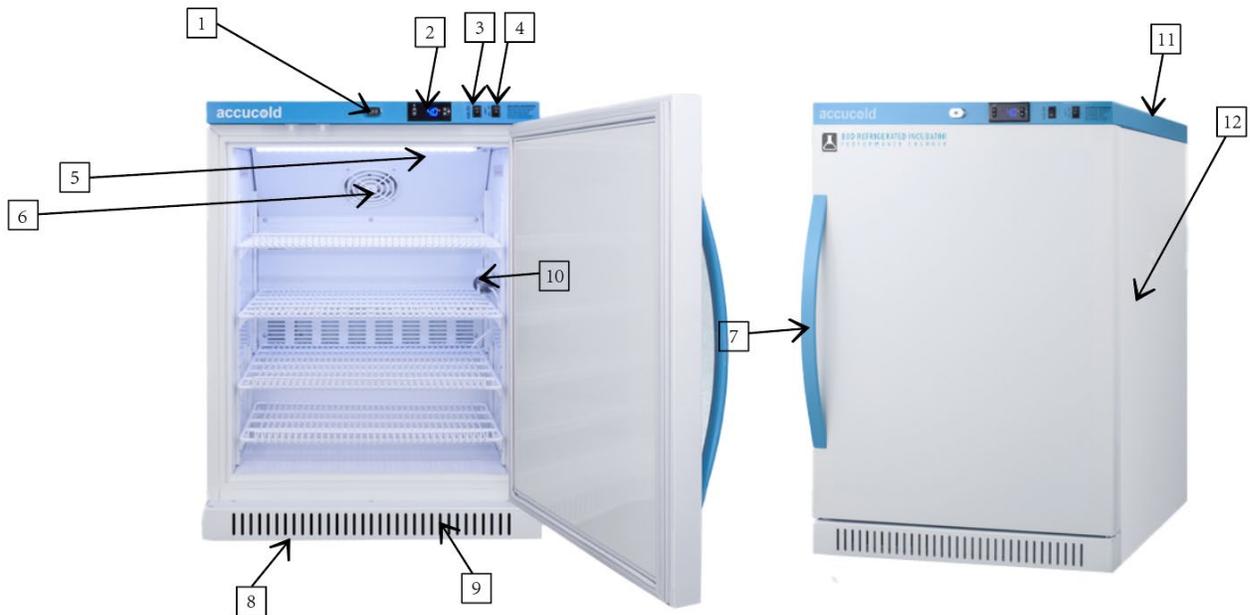

- | | |
|--|---|
| 1. Keyed Lock | 8. Leveling legs |
| 2. Microprocessor Control Panel | 9. Kickplate |
| 3. Reset Switch | 10. Temperature sensor in glycol bottle |
| 4. Light Switch | 11. Dry contacts (located in rear) |
| 5. Internal Light | 12. Access port – for added monitoring devices of the useable area inside the unit |
| 6. Internal Fan | |
| 7. Handle | |

Specification

Model	ABOD6	ABOD15	ABOD18
Working Temperature	0-38°C, RH 20-75%		
Operating Temperature	4-60°C (39.2 - 160 °F)		
Altitude	Up to 2000M		
Transport Temperature	-15°C – 50°C (5°F – 122°F)		
Storage Temperature	40-60°C / 104-140°F		
Capacity(L)	168	448	550
Rated Voltage	AC110V		
Mains Fluctuations	±10%		
Frequency	60Hz		
Rated Current	2.9A	4.7A	4.7A
Cooling Power	102W	200W	220W
Heating Power	300W	500W	500W
Shelf Load	20kg (44 lbs.) each		
Pollution Degree	2		
Product Dimensions MM (W*D*H)	595*611*840 23" x 24" x 33"	595*611*1820 23" x 24" x 72"	700*660*2020 28" x 26" x 80"
Net Weight	54.96 kg. 121.1 lbs.	90.5 kg 199.6 lbs.	123 kg. 269 lbs.
Refrigerant	R600a	R600a	R290
Filling Amount	1.34 oz	2.43 oz	3.0 oz
Vesicant	Cyclopentane		

OPERATION – QUICK SETUP GUIDE

This section provides an overview of how to operate your refrigerated incubator. **For more detailed instructions on adjusting control panel parameters, refer to page 10.** Once plugged in, the display will show the internal product temperature, and the compressor will begin running. The corresponding indicator light on the control panel will also turn on.

Sensors Overview:

The unit uses two sensors:

- **Product Sensor** (display sensor): Monitors the product temperature.
- **Air Sensor** (control sensor): Monitors the air temperature for controlling the refrigeration process.

For maximum performance, the air sensor is used to quickly respond to changes like product loading and door openings. You may need to adjust the controller to a lower temperature than the desired product temperature, as the air temperature needs to be lower to effectively cool the product.

The temperature displayed is the product temperature recorded by the product sensor inside a glycol-filled vial located in the unit. As the refrigerated incubator cools to the set point, the number on the display will decrease in temperature. It will take longer for the product vial to reach desired temperature compared to the air inside of the unit.

Power Outage


If power is cut off, wait for at least 5 minutes before plugging the unit in again to avoid damaging the compressor.

Control and Temperature Measurement Device (TMD)

Adjusting the Operating Temperature

The unit is preset to 20°C, adjust the temperature up or down till it reaches the desired product temperature. The displayed will reflect the current temperature setting.

To change the set point temperature, press the **SET** key. The set-point temperature will be displayed. Within 10 seconds, press the **UP** or **DOWN** arrow key to adjust the temperature. Press the **SET** key again to confirm the new setpoint.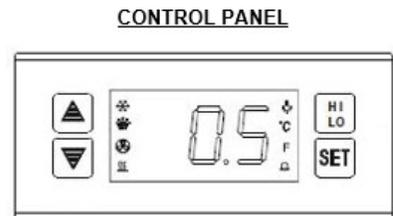

High/Low Temperature Alarm

- Low Temperature Alarm: If the air temperature in the refrigerated incubator falls below the low alarm set point, an alarm will sound. The front panel will display [LO] flashing alternately with the current temperature.
- High Temperature Alarm: If the air temperature exceeds the high alarm set point, an alarm will sound. The front panel will display [HI] flashing alternately with the current temperature.
- High Product Temperature Alarm: If the product temperature exceeds the high product alarm set point, the front panel will display [H2] flashing alternately with the current temperature.

Investigate the potential cause immediately if any of the alarms occur. The unit should be monitored to ensure that it returns to the correct temperature.

Adjust the alarms set points to meet your requirements in the parameter setup. (see page 13)

Power Failure Alarm

After plugging in the equipment, move the **Reset Switch** located at the front of the unit to the down position. When electrical power to the equipment is lost, an alarm will sound, and it will silence when the power is restored.

Note the alarm is battery-powered and has a limited time to alarm.

Min/Max Temperature Recording

Once connected to the power supply, your unit will record the maximum high and minimum low temperatures reached inside the cabinet. The display will show the last recorded **HI** and **LO** temperatures until the unit resets.

- To display the **HI** reading, press and release the **HI/LO** button.
- To display the **LO** reading, press and release the **HI/LO** button again.

To Reset the Min/Max Temperature Recording

Press and hold the **HI/LO** button for 6 seconds. To confirm the operation, The display will start blinking and the buzzer sounds. After 5 seconds, the display will revert to normal.

Changing the Temperature Display

The unit's default temperature is set to Celsius. If you prefer the temperature to display in Fahrenheit, follow these steps:

1. Enter Password Mode: Press and hold the **SET** button until the display flashes [PAS].
2. Enter the Password: Use the **UP** and **DOWN** arrows to enter the password (the factory default is 15).
3. Access the Parameter Settings: After entering the password, press the **SET** button about 25 times until [CF] appears on the display.
4. Change to Fahrenheit: Press the **DOWN** arrow until the display shows °F.
5. Save the Setting: Press **SET** and wait for 6 seconds. The new setting will be saved, and the unit will resume normal operation with the temperature displayed in Fahrenheit.

Interior Light

Your unit includes an LED light that illuminate automatically when the door is open. If you prefer the light to remain on when the door is closed, push the switch to the **ON** position.

Door Alarm

If the door is left open for more than 2 minutes, an alarm will sound. Close the door to stop the alarm.

Remote Alarm

The electronic controller is equipped with dry contacts for alarm forwarding to an external remote alarm system (via the terminal block on the back of the unit).

NOTE: The remote contacts are normally CLOSED and OPEN upon alarm. To change the output to normally OPEN and CLOSED on alarm see item **C12** of the "Table of Parameters and Factory Settings". Change the **C12** setting to **00** in lieu of **01**.

Probe Port

This unit is equipped with a 3/8" probe port and plug, which can accept additional monitoring devices. It is advisable to position the additional temperature monitoring device in the net usable area (inside ideally near the product sensor).


Summary of Alarm Types

ALARM TYPE	DISPLAY FLASH
Air temperature high alarm	HI delay alarm
Air temperature low alarm	LO delay alarm
Air temperature sensor short circuited	1H in time alarm
Air temperature sensor open circuited	1L in time alarm
Product temperature sensor short circuited	2H in time alarm
Air temperature sensor open circuited	2L in time alarm
Product temperature high temp alarm	H2 delay alarm
Product temperature low temp alarm	L2 delay alarm
Electric supply off alarm	No display – alarm in time
Door open delay alarm	Dr flash and delay alarm

Storage of Articles

- At maximum loading level, the content should not block the unit's air flow or be loaded above the load line.
- Leave space around the contents to allow a smooth flow of cold air inside the cabinet. To prevent freezing, be sure the articles do not touch the back of the interior cabinet.
- Store articles away from the inner fan.
- Shelves can be adjusted to allow proper airflow around the products being stored.

DETAILED TEMPERATURE CONTROLLER GUIDE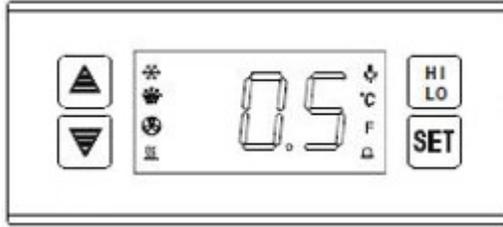

Front Panel Operation

Set Temperature (Compressor's Top Temperature) Adjustment:

1. Press the **SET** button, the set temperature is displayed, then press the **UP** or **DOWN** arrow to store and memorize. Press the **SET** button to exit the adjustment status and display the product temperature. If no other buttons are pressed within 6 seconds, the refrigerator product temperature will be displayed.
2. Set temperature adjustment range: Parameter E1~E2

Manually Start/Stop Defrosting:

Press **HI/LO** button, then press **SET** button and hold for 6 seconds to enter defrost status or to stop defrost.

Display The Sensor Temperature (C13):

Press the **DOWN** arrow, the product temperature (or cabinet temperature) will flash display. After 6 seconds, will resume normal display.

Refrigeration LED:

During refrigeration the cooling snowflake LED is on. When temperature is constant, the LED is off.

Defrost LED:

During defrosting (if activated), the defrost LED is on. During the delay after defrosting, the LED flashes.

Parameter Set-Up and Settings (For All Settings Other Than Temperature Setpoint):

Press **SET** button and hold for 6 seconds to enter the parameter setting (flash and display PAS). After entering the correct password (factory default is 15), press **SET** and the display will show E1, E2, ...~do3 PAS in sequence. Press the **UP** or **DOWN** arrow and the value of the parameter will be displayed and can be modified and stored. If no other buttons are pressed within 6 seconds, it will exit, and the new value will be stored.

NOTE: Only when in the inner parameter menu (display PAS) and the correct password is entered can the parameter value be adjusted. If the wrong password is entered, the parameter modification will be exited, although the set temperature adjustment is still active. If you forget the password, you will need to resume the factory default settings.

The following are examples of how to change common settings:

- A. **Enter Password** - Press and hold the **SET** button 6 seconds until it flashes [PAS] then enter the default password **15** using the **UP** and **DOWN** arrows. When **15** is reached press the **SET** button.
- B. **Change display to °F in lieu of °C** - After entering the password press the **SET** button about 25 times to reach the [CF] display, then press the up or down arrow until the control displays [°F] then press **SET**. Wait 6 seconds and it will be set and the unit will go back to normal operation.
- C. **High Air Alarm Settings** - After entering the password press the **SET** button until you reach the [C1] display, then press the **UP** or **DOWN** arrow until the control displays your desired setting then press **SET**. Wait 6 seconds and it will be set and the unit will go back to normal operation.
- D. **Low Air Alarm Settings** - After entering the password press the **SET** button until you reach the [C2] display, then press the **UP** or **DOWN** arrow until the control displays your desired setting then press **SET**. Wait 6 seconds and it will be set and the unit will go back to normal operation.
- E. **High Product Alarm Settings** - After entering the password press the **SET** button until you reach the [C3H] display, then press the **UP** or **DOWN** arrow until the control displays your desired setting then press **SET**. Wait 6 seconds and it will be set and the unit will go back to normal operation.
- F. **Low Product Alarm Settings** - After entering the password press the **SET** button until you reach the [C3L] display, then press the **UP** or **DOWN** arrow until the control displays your desired setting then press **SET**. Wait 6 seconds and it will be set and the unit will go back to normal operation.
- G. **Other Parameters in the "Factory Settings"**: Enter the Password and then press the **SET** button until the desired parameter is displayed, then press the **UP** or **DOWN** arrow until the desired setting is displayed then press the **SET** key again. Wait 6 seconds and it will be set and the unit will go back to normal operation.

Recording The Highest and Lowest Temperature

Once turned on and after the [C5] delay, the unit will start to record the highest and lowest temperature; the record will be refreshed at any time. The values will be kept in memory even after the power is turned off. Press the **HI/LO** button to see the highest temperature recorded. Press it again to see the lowest temperature. Holding the **HI/LO** button for 6 seconds will cause the display to flash and a buzzer to sound; the previous highest and lowest temperature readings will be cleared, and a new record will begin. (Highest and lowest temperature recording instructions: When parameter **F4=0**, the unit can record highest and lowest temperature at any time. When **F4=1,2,3**, during defrost and delay locking 20 minutes, highest and lowest temperatures are not recorded.

During air sensor failure, highest and lowest temperatures are not recorded. When set **C13=00**, the temperature in the liquid bottle is recorded; when set **C13=01**, the temperature of the air in the cabinet is recorded. The product temperature corresponds to the temperature of the liquid in the bottle.

Factory Default Resumption

Press and hold the **DOWN** arrow; at the same time, press the **UP** arrow and hold for 6 seconds. The display will flash and show [888]. At this time, all parameters will resume factory default values. After 6 seconds, normal operational mode will return.

To Check or Change the Password:

Enter the inner parameter menu (PAS display) and enter the correct password. After entering the inner parameter PAS, press the **UP** or **DOWN** arrow to display and change the password, then press **HI/LO** to confirm and store the new password.

Table of Parameters and Factory Settings

Parameter	Function	Set Range	Default	Parameter	Function	Set Range	Default
Set	Setpoint		20°C	C3H	Product Temp. High Temp. Alarm	C3L~80.0°C/176°F	70°C/158°F
PAS	Password	00~99	15	C3L	Product Temp. Low Temp. Alarm	-40°C/-40°F~ C3L	0.0°C/32.0°F
E1	Lowest set point limit	Set temp. ~ -4°C /39.2°F	4°C / 39.2°F	C4	Hysteresis Alarm	0.1~20.0°C / 0.2~36.0°F	0.1°C / 32.18°F
E2	Higher setpoint limit	60.0°C / 140.0°F ~set temp.	60.0°C / 140°F	C5	Start-up Temp. Alarm delay	00~99 min	60 min
E30	Temp. Hysteresis	1.5°C / 34.7°F	1.5°C / 34.7°F				
E31	Heating Temp. Hysteresis	0.1~20.0°C / 32.18~68.0°F	1.5°C / 34.7°F	C6	Temp. Alarm Delay	00~99 min	15 min
E40	Turning On delay time	00 ~10 min.	3 min.	C7	Power Off relay Alarm	00=do not alarm /01=alarm	01
E41	Comp. start delay time	00 ~10 min.	3 min.	C8	Alarm relay close after, muffle alarm relay switch	00=open / 01=close	00
E5	Offset on Air Temperature	-10.0~10.0°C / - 14.0~50.0°F	2.0°C / 35.6°F	C9	Restart time after buzzer mute	00=do not start / 01 ~30 min=restart time	10 min
E6	Offset on Product Temp.	-10.0~10.0°C / 14.0~50.0°F	0.0°C/ 32°F	C10	Comp. Force stop time	01~99 min	99 min
F0	Defrost Type	00=defrost by turning of comp.	00	C11	Comp. Force RunningTime	00=comp. Stop / 01-99min=starting time	0 min
F1	Max defrost duration	01~60 min.	20 min.	C12	Alarm Output Type	00=contact actuation when alarm / 01=contactdisconnect when alarm	01
F2	Defrost Interval (XXhr/time)	00~24 hr.	00	C13	Normal State temp.display type	00=liquid bottle temp / 01=cabinetair temp.	00
F4	Temp. displaying on screen when defrost	00= normal Temp; 01=latest Temp. 02=Fixed Temp. ; 03=showing "dEF" on screen	00	CF	TemperatureUnit	°C=Celsius / °F=Fahrenheit	°C
C0	Hot and cold conversion delay time	00-30min	02min	do1	Door open alarm	00=do not alarm /01=99min=delayed alarm	02 min
C1	High Temp Air Alarm	C2~80.0°C / 176.0°F	70.0°C / 158°F	do2	Comp. status whendoor open	0=stop / 01=original status	0 1
C2	Low Temp Air Alarm	-40°C / -40°F~C1	0.0°C / 32.0°F	do3	Heating control mode selection	00=Controlled by parameters 01 = PID intelligent control	00

FUNCTION DETAILS

Temperature Control: After turning on the unit and reaching the delay time (parameter E40), the compressor starts operating when the cabinet temperature is higher than (set temperature + hysteresis) and will be off when the cabinet temperature is lower than the set point temperature.

To protect the compressor, it cannot be re-started unless the time after the compressor stops is longer than the delay time (Parameter E41).

Defrost Control: After running for a defrost interval set time (parameter F2), the unit will enter defrost and the compressor stops. When the defrost duration time (parameter F1) ends, it will exit defrost status. After a defrost period ends, there will be 2 minutes of dripping time before the unit can enter refrigeration status.

- When defrost interval F2 is set to **00**, the automatic defrost by turning off compressor will be cancelled.
- When setting parameter F4 is set to **0**, the temperature will be displayed normally during defrost.
- When setting parameter F4 is set to **1**, the cabinet temperature is locked during defrost, and the last value before defrost is displayed. When defrost ends, normal display will resume after temperature display 20 minutes delay (or cabinet temperature lower than the set temperature). The defrost LED flashes during delay.
- When setting parameter F4 is set to **2**, the set temperature will be displayed during defrost.
- When defrost ends, normal display will be resumed after 20 minutes delay (or cabinet temperature lower than the set temperature). The defrost LED flashes during delay.
- When setting parameter F4 is set to **3**, [dEF] will be displayed during defrost. When defrost ends, normal display will be resumed after 20 minutes delay display [dEF] (or cabinet temperature lower than the set temperature). The defrost LED flashes during delay.

Alarm Control: After turning on for the first time, the unit will need to reach the [C5] delay time setting, then the high/low temperature alarm function can be triggered (C1, C2, C3). After passing [C5] delay, when the cabinet temperature is abnormal (for example higher than the temperature alarm [C1] or lower than temperature alarm [C2]) and duration more than alarm delay time [C6], the unit will enter alarm status, alarm start. When high temperature alarm alternate display H1 and cabinet air temperature, the compressor starts to run. When low temperature alarm, will alternate display **LO** and cabinet temperature, the compressor stops. When cabinet temperature is higher than the (low temperature alarm value [C2] + alarm hysteresis [C4]), the low temperature alarm ends. When cabinet temperature is lower than (high temperature alarm value [C1] – alarm hysteresis [C4]), the high temperature alarm ends.

When the product sensor temperature is greater than or equal to the product temperature high temperature alarm value [C3] and duration more than alarm delay time [C6], unit enters alarm status, and starts the alarm, alternate display [H2] and cabinet temperature. When the product sensor temperature is less than or equal to the (high temperature alarm value [C3] – alarm hysteresis [C4]), it will end the high temperature alarm.

When the electric supply is off, flash and display [EEL] and alarm. When setting **do1** is set at **00**, unit will not alarm when door is open. When setting **do1** is set at **0**, once unit reaches the delay time, it will flash and display [dr alarm], press any button to cancel.

If setting [C7] is set to **0**, then the units relay does not alarm when power is off.

If setting [C8] is set to **0**, then after alarm is canceled, the alarm relay does not close. If [C9] is set to **0**, then after the alarm cancel button is pushed, the buzzer does not restart.

If setting other numbers, then after reaching delay time, buzzer sounds one more time. (Under the condition of alarm not terminating)

Abnormal Work Mode: When air sensor is short-circuited or high temperature over limit ($> 66^{\circ}\text{C}/151^{\circ}\text{F}$), [1H] is displayed; when air sensor is open-circuited or low temperature over limit ($< -40^{\circ}\text{C}/^{\circ}\text{F}$) [1L] is displayed. Compressor will enter the force running mode, according to C10, C11 set parameters running in sequence.


When product sensor is short-circuited, open-circuited or over limit ($> 66^{\circ}\text{C}/151^{\circ}\text{F}$), alternate display 2H and air temperature, the product temperature sensor open-circuited or low temperature over limit ($< -40^{\circ}\text{C}/^{\circ}\text{F}$), alternate display [2L] and air temperature.

Instrument Running Alarm Indication List

Alarm type	Compressor running	Display flash
Refrigerated incubator high temp. alarm	Compressor on	H1 delay alarm
Refrigerated incubator low temp. alarm	Compressor off	L0 delay alarm
Air temp. sensor short-circuited	Press C10, C11 to run the comp.	1H in time alarm
Air temp. sensor open-circuited	Press C10, C11 to run the comp.	1L in time alarm
Product temp. sensor short-circuited		2H in time alarm
Product temp. sensor open-circuited		2L in time alarm
Product temperature high temp.	Limit alarm	H2 delay alarm
Product temperature low temp.	Limit alarm	L2 delay alarm
Electric supply off alarm		EEL flash and alarm in time
Door open delay alarm		Dr flash and delay alarm

When the door is open, the light switch can turn the light ON and OFF.

Circuit Diagram

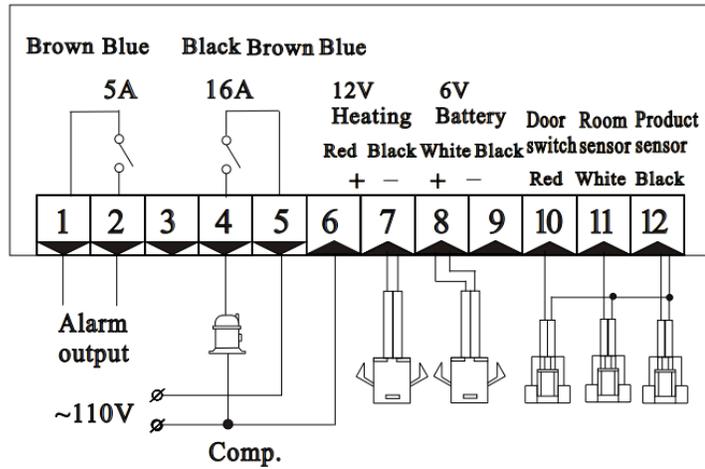- When the door is closed, normal close contact
- When the door is open, contact is disconnected

Notes for Installation:

1. Sensor leads must be kept separately from main voltage wires in order to avoid high frequency noise induced. Separate the power supply of the loads from the power supply of the controller.
2. When installing, the sensor must be placed with the head upward and the wire downward.
3. The temperature controller must not be installed in an area where water drips.

CLEANING AND MAINTENANCE

Before cleaning, it is essential that you unplug the unit from the power line and transfer the contents where they can be stored and monitored at the correct temperatures.

- Wash the inner compartment with warm water and neutral detergent. DO NOT allow the control panel, cables or plug to get wet. NEVER use corrosive detergents, wire brushes, or abrasive scourers to clean your refrigerated incubator. NEVER use metal or sharp implements to remove debris.
- Dry all surfaces thoroughly.
- To ensure trouble-free operation, the condenser should be cleaned every three months where appropriate using a vacuum hose. The condenser is located at the back of the cabinet. In exceptionally dusty locations, the condenser should be cleaned more often.
- When transporting the unit, try to keep it vertical with angle of inclination between sides of the refrigerated incubator and the horizontal plane not less than 50°. Otherwise the compressor may be affected, disrupting normal operation of the unit.

TROUBLESHOOTING

You can solve many common problems easily, saving you the cost of a possible service call. Try the suggestions below to see if you can solve the problem before calling the servicer.

Problems	Possible causes	Remedy
<i>The unit does not operate</i>	<i>Bad connection of plug or burnt out fuse</i>	<i>Unplug the power cord and re-connect the plug to power supply or replace with a new fuse</i>
<i>Abnormal operation of the compressor or there is a buzzing sound</i>	<i>The power voltage is out of rated range</i>	<i>Disconnect the power supply immediately and reconnect with it after normal voltage. It is necessary to have a voltage stabilizer in case of poor power supply.</i>
<i>Compressor operates for a long time and no frost on the surface of the evaporator</i>	<i>Refrigeration system is at fault (leakage or blocked)</i>	<i>Call for service</i>
<i>There is frost or ice on the walls of the inner cabinet and internal temperature is too low, the compressor never stops running.</i>	<i>Thermostat does not work</i>	<i>Call for service</i>
	<i>The temperature setting of the thermostat is too low.</i>	<i>Adjust to a higher temperature</i>
<i>The internal temperature is too high, and the compressor never stops operating.</i>	<i>Bad heat dissipation and ventilation of condenser</i>	<i>Improve ventilation</i>
	<i>Too many warm items were put in at one time</i>	<i>Remove some goods so air can circulate</i>
	<i>Door is being opened too frequently during initial cool down</i>	<i>Permit the unit to cool down adequately, prevent product access during this</i>
<i>Too noisy</i>	<i>The unit is not level</i>	<i>Adjust the adjustable legs</i>
	<i>The fastener of the unit is loose</i>	<i>Tighten the loose fastener</i>
	<i>Pipe near the compressor are touching</i>	<i>Carefully separate the touching pipes</i>
<i>The side of unit is hot</i>	<i>The condenser in the side wall gives out heat as part of normal operation</i>	<i>Nothing to worry about.</i>
<i>Sometimes a light sound of water flowing will be heard</i>	<i>Refrigerant flowing inside the pipe.</i>	<i>Nothing to worry about.</i>
<i>There may be condensation on the glass door</i>	<i>High ambient temperature or humid conditions</i>	<i>Dry with a cloth</i>
<i>The display shows "So" and compressor does not run</i>	<i>The temperature sensor is open circuit</i>	<i>Call for service</i>
<i>The display shows "SC" and compressor does not run</i>	<i>The temperature sensor is short circuited.</i>	<i>Call for service</i>

<i>The display shows "HH" and compressor does not run</i>	<i>The ambient temperature is above 40°C/104°F</i>	<i>Stop using immediately</i>
<i>The display shows "LL" and compressor does not run</i>	<i>The ambient temperature is below -40°C/-40°F.</i>	<i>Stop using immediately</i>

If you continue to experience issues, contact Accucold's technical support department at **1-888-4-MEDLAB**.

TECHNICAL INFORMATION

Model Numbers	Gross Capacity Cu. Ft.	Net Useable Interior Capacity Cu. Ft.
ABOD6	6.00	4.40
ABOD15	15.00	12.10
ABOD18	18.00	17.40

LIMITED WARRANTY

TWO-YEAR LIMITED WARRANTY

Within the 48 contiguous United States, for two years from the date of purchase, when this unit is operated and maintained according to instructions attached to or furnished with the product, warrantor will pay for factory- specified parts and repair labor to correct defects in materials or workmanship. Service must be provided by a designated service company. Outside the 48 states, all parts are warranted for two years from manufacturing defects. Plastic parts, shelves and cabinets are warranted to be manufactured to commercially acceptable standards and are not covered from damage during handling or breakage.

FIVE-YEAR COMPRESSOR WARRANTY

1. The compressor is covered for five years.
2. Replacement does not include labor.

ITEMS WARRANTOR WILL NOT PAY FOR:

1. Service calls to correct the installation of your equipment, to instruct you how to use your equipment, to replace or repair fuses or to correct wiring or plumbing.
2. Service calls to repair or replace unit light bulbs or broken shelves. Consumable parts (such as filters) are excluded from warranty coverage.
3. Damage resulting from accident, alteration, misuse, abuse, fire, flood, acts of God, improper installation, installation not in accordance with electrical or plumbing codes, or use of products not approved by warrantor.
4. Replacement parts or repair labor costs for units operated outside the United States.
5. Repairs to parts or systems resulting from unauthorized modifications made to the unit.
6. The removal and reinstallation of your unit if it is installed in an inaccessible location or is not installed in accordance with published installation instructions.

DISCLAIMER OF IMPLIED WARRANTIES; LIMITATION OF REMEDIES CUSTOMER'S SOLE AND EXCLUSIVE REMEDY UNDER THIS LIMITED WARRANTY SHALL BE PRODUCT REPAIR AS PROVIDED HEREIN. IMPLIED WARRANTIES, INCLUDING WARRANTIES OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE, ARE LIMITED TO ONE YEAR. WARRANTOR SHALL NOT BE LIABLE FOR INCIDENTAL OR CONSEQUENTIAL DAMAGES. SOME STATES DO NOT ALLOW THE EXCLUSION OR LIMITATION OF INCIDENTAL OR CONSEQUENTIAL DAMAGES, OR LIMITATIONS ON THE DURATION OF IMPLIED WARRANTIES OF MERCHANTABILITY OR FITNESS, SO THESE EXCLUSIONS OR LIMITATIONS MAY NOT APPLY TO YOU. THIS WARRANTY GIVES YOU SPECIFIC LEGAL RIGHTS AND YOU MAY ALSO HAVE OTHER RIGHTS, WHICH VARY FROM STATE TO STATE.

WARNING: This product can expose you to chemicals including Nickel (Metallic) which is known to the State of California to cause cancer.

For more information go to www.P65Warnings.ca.gov

Note: Nickel is a component in all stainless steel and some other metallic compositions.

Accucold Division of Felix Storch, Inc.
An ISO 9001:2015 registered company770
Garrison Avenue
Bronx, NY 10474

accucold[®]
For parts and accessory ordering,
troubleshooting and helpful hints, visit:
www.accucold.com/support

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